

Manual for RoR cleaning set #206

12-jan-2025

- 1)First blow away as much loose dust as possible with the GIOTTO blow balloon.
- 2)Take the bottle of RoR-1 lens cleaner and first moisten the tip of a cleaning stick, then twist a piece of cotton wool on the end of the stick (without touching it with your fingers. This is best done by turning the stick in your elbow on the fabric of your (clean) sleeve.
- 3)Now moisten the cotton wool tip with a little RoR.
- 4)Place the objective upright on the table and turn the cotton wool tip around on the front lens without applying force. To dry; repeat 2) but now without RoR.
- 5)Check for cleanliness with an eyepiece as a magnifying glass (held backwards).
- 6)Repeat the procedure in point 3 (always use a new piece of cotton wool!
- 7)Always check with the eyepiece, and repeat these procedures until the front lens is spotlessly clean.

IN GENERAL

- Never use too much force when cleaning.
- Always use a clean piece of cotton wool.
- Use a small amount of cleaning agent
- In case of stubborn dirt, repeat the procedure after a while, if -necessary moisten it first with RoR and wait a while.

RoR lens cleaner is an extremely safe product for cleaning of:

Objectives-Eyepieces-CCDs-Laserlenses-Flatscreens-Glasses-Monitor-Mobile phones-Scanners-Telescopes-Binoculars-Filters-Aiming sights-Safety glasses-Projection lenses etc.etc.